

Preservation  
above-zero



MPR-S163

Temperature range

**+2°C to 14°C**

**Compact and large capacity**

**New cycle defrosting system**

### Medical Refrigerators for reliable storage of pharmaceutical and medical supplies

- Precise temperature control
- Superior cooling performance
- Forced air circulation
- Double-glazing glass door
- Slim, and space saving design
- Useful alarm functions

# MPR-S163

## Features

### Improved Energy Saving

PHCbi's new cycle defrosting system achieved 15 %\* energy saving compared to conventional models. This system observes the defrosting condition after every cycle, and prevents temperature rises that are caused by defrosting.

\*The data is a measured example (set temp. 5°C, AT 23°C, 230 V 50 Hz, no load). Actual power consumption will vary depending on the set and ambient conditions, loading and local voltage.

### User Friendly Design

#### 1. New Control Panel

- Newly designed for improved operability, visibility, and calibration
- Digital setup for easy and secure operation
- Temperature alarm settings made on the control panel



#### 2. LED Interior Light

The LED interior light automatically turns on/off in conjunction with the door opening/closing. It can also be controlled from the control panel.



#### 3. Double-Paneled Glass with Heat-Reflective Film

The refrigerator doors constructed of double-paneled glass with heat-reflective film, allow easy loading and unloading of medical supplies while preventing ultraviolet rays, which may damage stored items, from entering the unit. Panasonic's unique heat-reflective film blocks the passage of radiant heat rays through the glass panels and keeps the inside temperature from being adversely affected by excessive amounts of heat.

## Effective Temperature Control

### 1. Microprocessor Temperature Control

A thermistor sensor monitors temperature inside the chamber, and microprocessor temperature control ensures that the set temperature is maintained. Even if the door is opened and closed frequently, the circulation fan ensures rapid temperature adjustment to provide a highly reliable, stable preservation environment that is not affected by ambient temperature.

### 2. Fan-Forced Air Circulation

The temperature stays even throughout the inside of the refrigerator with the fan-forced air circulation system. No matter how the load is distributed, every corner of the unit is immediately cooled with no noticeable variation in temperature apparent inside the cabinet.

## Enhanced Alarm and Safety Functions

### 1. Door Open Alarm

When the door is opened, the Door Open indicator lamp lights automatically. After approx. 2 minutes, a buzzer alarm will sound if the door has not been closed (settable from 0 to 15 minutes). If the door remains open even after the buzzer is stopped, the buzzer rings again after 30 minutes (settable from 0 to 60 minutes in increments of 10 minutes).

### 2. Abnormal Temperature Alarm

If the inside temperature changes  $\pm 2^{\circ}\text{C}$  to  $\pm 14^{\circ}\text{C}$  from the set temperature, the digital temperature display flashes, and a buzzer sounds an alarm after 15 minutes (settable from 0 to 15 minutes).

## Specifications

	Model No.
220V-240V, 50Hz	MPR-S163-PE
220V, 60Hz	MPR-S163-PK MPR-S163-PR MPR-S163-PT
110/115V, 60Hz	—
230/240V, 50Hz	—
220V, 50Hz/60Hz	—
<b>External dimensions (W x D x H)*1</b>	800 x 465 x 1090 (mm) 31.5 x 18.3 x 42.9 (inch)
<b>Internal dimensions (W x D x H)</b>	720 x 300 x 725 (mm) 28.3 x 11.8 x 28.5 (inch)
<b>Effective capacity</b>	158 liters (5.6 cu.ft.)
<b>Net weight</b>	71 kg (156 lbs.)
<b>External cabinet</b>	Galvanised steel with baked-on finish
<b>Internal cabinet</b>	Stainless steel
<b>Insulation</b>	Polyurethane foam
<b>Doors</b>	Sliding glass doors, double-glazing glass with heat-reflective film
<b>Shelves</b>	Hard steel wire
<b>Access port</b>	30mm $\phi$ on back wall
<b>Lighting / Casters</b>	LED/2 casters
<b>Compressor</b>	Hermetic type, 90 W
<b>Refrigerant</b>	HFC
<b>Evaporator</b>	Fin & Tube, Forced-air circulation
<b>Condenser</b>	Wire & Tube
<b>Defrosting</b>	Cyclical defrosting & evaporator temp. detection system
<b>Defrosting heaters</b>	87 W
<b>Temperature control</b>	2°C to 14°C
<b>Temperature display</b>	Digital (1°C increments)
<b>High/Low temperature alarm system</b>	$\pm 2^{\circ}\text{C}$ to $\pm 14^{\circ}\text{C}$ from temperature setting value
<b>Door ajar alarm</b>	Buzzer / door ajar lamp

**Options** Temperature recorder: MTR-0621LH-PE (recorder fixing: MPR-S30-PW)  
MTR-G04C-PE (recorder fixing: MPR-S7-PW)

Battery mounting box: MPR-48B1-PW

Interface board (For data acquisition system MTR-5000 users only)

MTR-L03-PW / MTR-480-PW


**Intended use:** This equipment is designed for storage of pharmaceuticals, samples and reagents.

**Caution:** PHC Corporation guarantees this product under certain warranty conditions. However, please note that PHC Corporation shall not be responsible for any loss or damage to the contents of the product.

- Appearance and specifications are subject to change without notice.
- Water condensation may sometimes appear on the glass doors, but this is normal and not a malfunction.

\*1 External dimensions of main cabinet only - see dimension drawings showing handles and other external projections.

## Dimensions Unit: mm (inch)

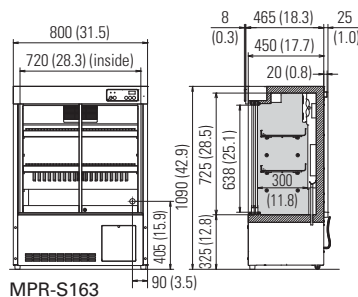

Preservation Equipment, Experimental Environment Equipment, Dispensary Equipment, Culturing Equipment and Drying & Sterilising Equipment for General Laboratory use  
The management of the design, development, production and servicing of the above.  
PHC Corporation, Biomedical Division  
1-1-1 Sakada, Oizumi-machi, Ora-gun, Gunma 370-0596, Japan



PHC Corporation, Biomedical Division is certified for:  
**Environmental management system: ISO14001**

## DISTRIBUTED BY:

 **PT. SIGMA BIMED**

**Indonesia Sole Distributor**

Jl. Agung Utara Raya Blok A 36D/42 Sunter Agung Jakarta Utara 14350 Indonesia

Teip. (62-21) 64716568, 70240088, 70240099, 87960123, 87960456

Fax. (62-21) 64716566, 87961123, 87961456

Email : sigmabimed@indo.net.id, sales@sigmabimed.com

Website : www.sigmabimed.com

Kemenkes RI AKL 20209914675

**phcbi**

**PHC Corporation**

1-1-1 Sakata Oizumi-Machi, Ora-Gun, Gunma 370-0596, Japan

© PHC Corporation